## Buffered NaCl-Peptone Solution modified pH 7.0 + NT

Diluent with neutralizers for detection and enumeration of microorganisms, according to USP/EP/JP and ISO 21149.

### DESCRIPTION

Buffered NaCl-Peptone Solution pH 7.0 modified + NT is used for dissolving, suspending and diluting test samples.

Its composition is based on the recommendations in the Harmonized Pharmacopoeias for the microbiological examination of non-sterile products.

Conforms to ISO 21149 for the detection of viable microorganisms in cosmetics, this formula enriched with neutralizers allows the preparation of samples which have an antimicrobial effect or which are not soluble in water.

### TYPICAL FORMULA

	(g/l)
Potassium Dihydrogen Phosphate	3.6
Disodium Hydrogen Phosphate Dihydrate	7.2
Sodium Chloride	4.3
Peptone	1.0
Lecithin	3.0
Histidine	1.0
Polysorbate 80	30 ml
Sodium Thiosulfate	5.0
Final pH 7.0 ± 0.2 at 25°C	

### METHOD PRINCIPLE

Sodium phosphate and potassium phosphate are the buffering agents. Sodium chloride maintains the osmotic balance. A low peptone content provides nitrogen, vitamins, minerals and amino acids to ensure the viability of the organisms. Lecithin neutralizes quaternary ammonium compounds. Histidine inactivates aldehydes. Polysorbate (Tween 80) is a source of fatty acids and is effective against phenolic compounds and mercurial derivatives. Sodium thiosulfate neutralizes halogen compounds.

### TEST PROCEDURE

Dissolve or dilute the product to be examined (usually a 1 in 10 dilution is prepared) in Buffered NaCl-Peptone Solution modified pH 7.0 + NT. Use the suspensions within 2 hours or within 24 h if stored at 2-8°C.

### INTERPRETING RESULTS

No marked increase or decrease in original colony forming unit (CFU) count.

### APPEARANCE

Clear, colorless.

### STORAGE

Store tubes and bottles at 10-25°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

### SHELF LIFE

2 years.

**QUALITY CONTROL**

The medium is inoculated with the microbial strains indicated in the QC table.

Inoculum for diluent:  $10^3$ - $10^4$  CFU.

Incubation conditions:  $22.5 \pm 2.5^\circ\text{C}$  for 2 hours.

**QC Table.**

Microorganism		Growth on TSA
<i>Staphylococcus aureus</i>	ATCC® 6538	$\pm 30\%$ colonies of original count
<i>Pseudomonas aeruginosa</i>	ATCC® 9027	$\pm 30\%$ colonies of original count
<i>Escherichia coli</i>	ATCC® 8739	$\pm 30\%$ colonies of original count
<i>Salmonella Typhimurium</i>	ATCC® 14028	$\pm 30\%$ colonies of original count
<i>Bacillus subtilis</i>	ATCC® 6633	$\pm 30\%$ colonies of original count
<i>Candida albicans</i>	ATCC® 10231	$\pm 30\%$ colonies of original count

**WARNING AND PRECAUTIONS**

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

**DISPOSAL OF WASTE**

Disposal of waste must be carried out according to national and local regulations in force.

**BIBLIOGRAPHY**

1. EN ISO 11133:2014. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media.
2. European Pharmacopoeia 6.5 (2009) 2.6.13. Microbiological examination of non-sterile products: Test for specified microorganisms.
3. United States Pharmacopoeia 32 NF 27 (2009) <62> Microbiological examination of non-sterile products: Test for specified microorganisms.
4. Japanese Pharmacopoeia 4.05 (2008) Microbiological examination of non-sterile products: Test for specified microorganisms.
5. ISO 21149:2006 Cosmetics - Enumeration and detection of aerobic mesophilic bacteria.

**PRESENTATION**

	Contents	Ref.
Buffered NaCl-Peptone Solution pH 7.0 modified + NT	100 x 9 ml tubes	26470
Buffered NaCl-Peptone Solution pH 7.0 modified + NT	25 x 90 ml bottles	452650

**TABLE OF SYMBOLS**

<b>LOT</b> Batch code	 Keep away from sunlight	 Manufacturer	 Use by	 Fragile, handle with care
<b>REF</b> Catalogue number	 Temperature limitation	 Contains sufficient for <n> tests	 Caution, consult Instruction For Use	 Do not reuse

**LIOFILCHEM® s.r.l.**

Via Scozia zona ind.le, 64026 Roseto degli Abruzzi (Te) Italy  
Tel. +39 0858930745 Fax +39 0858930330 www.liofilchem.net liofilchem@liofilchem.net

## Buffered NaCl-Peptone Solution modified pH 7.0 + NT

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