# Baird Parker Agar Base + RPF Supplement

Selective medium for the isolation of coagulase-positive staphylococci in food, according to ISO 6888-2.

## DESCRIPTION

Baird Parker Agar Base + RPF Supplement is a kit used for preparing a selective and differential medium intended for the isolation and enumeration of coagulase-positive staphylococci from food and other materials, according to ISO 6888-2.

## KIT CONTENTS

Each kit contains:

- 6 bottles of Baird Parker Agar Base
- 6 vials of lyophilized RPF Supplement
- 1 instruction sheet

## PRINCIPLE OF THE METHOD

Pancreatic digest of casein, yeast extract and meat extract provide carbon, nitrogen, amino acids, vitamins and minerals for organisms growth. Sodium pyruvate and glycine stimulate growth of staphylococci. Agar is the solidifying agent. Lithium chloride and potassium tellurite make the medium selective for staphylococci by suppressing the accompanying flora. Fibrinogen and rabbit plasma allow the coagulase reaction. Trypsin inhibitor is added to prevent fibrinolysis.

## COMPOSITION

Baird Parker Agar Base	Content / liter of medium	RPF Supplement	Content / vial
Pancreatic Digest of Casein	10.0 g	Potassium Tellurite	2.5 mg
Yeast Extract	1.0 g	Bovine Fibrinogen	0.5 g
Meat Extract	5.0 g	Rabbit Plasma – EDTA	2.5 ml
Sodium Pyruvate	10.0 g	Trypsin Inhibitor	2.5 mg
L-Glycine	12.0 g		
Lithium Chloride	5.0 g		
Agar	17.0 g		
Final pH 7.2 ± 0.2			

## PROCEDURE FOR USE

Melt the content of 1 bottle of Baird Parker Agar Base in a boiling water bath at 100°C (loosing the caps partially unscrewed) until completely dissolved. Cool down to 48 ± 1°C. Aseptically, add the content of 1 vial of RPF Supplement reconstituted with 10 ml of sterile distilled water. Mix well avoiding foam formation.

## TECHNIQUE AND INTERPRETATION OF THE RESULTS

ISO 6888-2 recommends to inoculate the medium by using the Pour Plate Method. Incubate plates at 37 ± 1°C for 24 ± 2 hours. If necessary, re-incubate for 24 ± 2 hours. Staphylococci form black or grey or even white, small colonies surrounded by halo of precipitation, indicating coagulase activity.

## QUALITY CONTROL

1. Control of the appearance: amber medium, slightly opalescent. Freeze-dried pinkish-grey in colour.
2. Microbiological control.

Prepare the medium as indicated in the procedure for use. Inoculate the plates with the strains indicated below and incubate at 37 ± 1°C for 24-48 h.

Control strains	WDCM	Growth	Specification
<i>Staphylococcus aureus</i>	WDCM 00034	Good	Black or grey colonies with opacity halo
<i>Escherichia coli</i>	WDCM 00013	Inhibited	---
<i>Staphylococcus saprophyticus</i>	WDCM 00159	Good	Black or grey colonies without opacity halo
<i>Staphylococcus epidermidis</i>	WDCM 00036	Good	Black or grey colonies without opacity halo

## PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

## STORAGE

2-8°C in its original packaging. Keep away from sources of heat and avoid excessive changes of temperature. Use until the expiry date indicated on the label. Eliminate without using if there are signs of deterioration. Once reconstituted, the product can be stored for a maximum duration of 30 days at -20°C, shielded from light.

## SHELF LIFE

2 years.

## REFERENCES

- ISO 6888-2 (1999) Microbiology of food and animal feeding stuffs – Horizontal method for the enumeration of coagulase-positive staphylococci (*Staphylococcus aureus* and other species) – Part 2: Technique using rabbit plasma fibrinogen agar medium.
- Baird Parker A.C. (1962) An improved diagnostic and selective medium for isolating coagulase-positive staphylococci. J. Appl. Bact. 25:12-19.

## PRESENTATION

Product	Ref.	Contents
Baird Parker Agar Base + RPF Supplement	420010	<ul style="list-style-type: none"> <li>• 6 x 90 ml bottles of Baird Parker Agar Base</li> <li>• 6 vials of RPF Supplement</li> </ul>

One vial of RPF Supplement is sufficient to prepare 100 ml of medium.

## TABLE OF SYMBOLS

<b>LOT</b>	Batch code	 Do not reuse	 Manufacturer	 Contains sufficient for <n> tests	 Temperature limitation
<b>REF</b>	Catalogue number	 Fragile, handle with care	 Use by	 Caution, consult accompanying documents	

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