# Eugon LT 100 Supplement

Liquid supplement for preparing Eugon Broth LT 100.

## DESCRIPTION

Eugon LT 100 Supplement is a liquid supplement containing Octoxynol 9 (Triton X-100) and Polysorbate 80 (Tween 80). Supplement is added along with Lecithin Supplement (ref. 80007) to Eugon Broth (Ref. 610341). The supplemented medium is used for the enrichment and detection of aerobic bacteria in cosmetic products.

## KIT CONTENTS

Each kit contains:

- 4 bottles of liquid Eugon LT 100 Supplement
- instruction sheet

## PRINCIPLE OF THE METHOD

Triton X-100 is an excellent detergent that enhances the release of microorganisms from the slack matrix of the cosmetic emulsion. Lecithin and Tween 80 are neutralizers that inactivate preservatives like quaternary ammonium compounds, phenol and aldehydes derivatives

## COMPOSITION

	Content / bottle	Content / liter of medium
Octoxynol 9 (Triton X-100)	6.7 g	1.0 g
Polysorbate 80 (Tween 80)	33.3 g	5.0 g

## PROCEDURE FOR USE

1. Aseptically add 3 ml Eugon LT 100 Supplement and 5 ml of Lecithin Supplement to 500 ml of Eugon Broth autoclaved and cooled at 45-50°C
2. Mix with care and pour into final containers.

## TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical sheet of the medium being prepared.

## QUALITY CONTROL

1. Visual inspection: dense and oily, amber liquid.
2. Microbiological control.  
Prepare the medium per label directions with the addition of supplements. Inoculate the broth with the microbial strains indicated below and incubate at  $32.5 \pm 2.5^\circ\text{C}$  for 24-72 h.

### Control strains

*Staphylococcus aureus* ATCC® 25923  
*Candida albicans* ATCC® 10231

### Growth

Good  
Good

## WARNING AND PRECAUTIONS

The product contains hazardous substances and is classified as dangerous. Consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

## STORAGE AND TRANSPORT CONDITIONS

2-8°C away from light, until the expiry date on the label. However, our stability studies have shown that the storage or transport at 18-25°C for 4 days, or at 35-39°C for 48 hours, do not alter in any way the performance of the product. Eliminate if signs of deterioration or contamination are evident.

## REFERENCES

- Guisno, R., I.W. Gibby & M.J. Foter (1946) A neutralizing medium for evaluation of the germicidal potency of the quaternary ammonium salts. Amer J Pharm 118:320-323.
- Pelczar and Vera (1949) Milk Plant Monthly 38:30.
- ISO 21149 Standard (2006) Cosmetics – Microbiology – Enumeration and detection of aerobic mesophilic bacteria.
- ISO 21150 Standard (2006) Cosmetics – Microbiology – Detection of *Escherichia coli*.

## PRESENTATION

Product	Ref.	Content
Eugon LT 100 Supplement	80048	4 x 40 ml bottles

One bottle is sufficient to prepare more than 6 liters of medium.

## TABLE OF SYMBOLS

 <b>LOT</b> Batch code	 Do not reuse	 Manufacturer	 Contains sufficient for <n> tests	 Temperature limitation
 <b>REF</b> Catalogue number	 Fragile, handle with care	 Use by	 Caution, consult accompanying documents	

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LIOFILCHEM® S.r.l.