MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement

Selective supplement for the detection of *Listeria spp.* from food.

DESCRIPTION

MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement is a selective supplement for the detection of *Listeria spp.* from food when used with BUFFERED LISTERIA ENRICHMENT BROTH code 610324 or 620324.

KIT CONTENTS

Each kit contains:

- 10 bottles of **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement**
- 1 Instruction sheet

PRINCIPLE OF THE METHOD

The composition in antibiotics guarantees the selectivity: nalidixic acid inhibits Gram-negative micro-organisms and acriflavine inhibits some Gram-negative and Gram-positive micro-organisms; Amphotericin B avoids fungi growth.

COMPOSITION

| MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement | | |
|---|------------------|-----------------------|
| | Content / bottle | Content / L of medium |
| Nalidixic acid | 20.0 mg | 40.0 mg |
| Amphotericin B | 50.0 mg | 100.0 mg |
| Acriflavine hydrochloride | 7.5 mg | 15.0 mg |

PROCEDURE FOR USE

Reconstitute the contents of a bottle of **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement** aseptically with 5 ml of sterile distilled water. Shake until completely dissolved, avoiding foam formation.

1. Add the entire contents of a bottle (5 ml) aseptically to 500 ml of BUFFERED LISTERIA ENRICHMENT BROTH code 610324 or 620324, autoclaved and cooled to 45-50°C.
2. Mix with care.
3. Distribute onto sterile containers.

TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical documentation of BUFFERED LISTERIA ENRICHMENT BROTH code 610324 or 620324.

QUALITY CONTROL

1. Control of the appearance: a white freeze-dried product.
2. Microbiological control.

Prepare the tubes using as medium BUFFERED LISTERIA ENRICHMENT BROTH code 610324 or 620324, enriched with **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement** (1 bottle in 500 ml of medium). Inoculate tubes with the strains indicated in the microbiological control table. Incubation conditions: 18-48 h at 30 ± 2 °C.

Microbiological control

| Micro-organisms | | Growth |
|-------------------------------|------------|-----------|
| <i>Listeria monocytogenes</i> | ATCC 35152 | good |
| <i>Enterococcus faecalis</i> | ATCC 29212 | inhibited |

PRECAUTIONS

The product **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement** is classifiable as hazardous under current legislation; it is recommended that the Safety Data Sheet be consulted on its use.

MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement must be used in the laboratory by properly trained personnel, using approved asepsis and safety methods for handling pathogenic agents.

STORAGE

Store **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement** at 2-8°C in its original packaging. Keep away from sources of heat and avoid excessive changes of temperature. In such conditions **MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement** maintains its validity until the expiry date indicated on the label. Eliminate without using if there are signs of deterioration.

REFERENCES

1. Lovett J., Francis D.W. and Hunt J.M. (1987) *J.Food Prot.* 50. 188-192.
2. Curtis G.D.W., Nichols W.W. And Falla T.J. (1989). *Lett Appl. Microb.* 8. 169-172.

PRESENTATION

| Product | REF | Σ |
|--|--------------|-------------------|
| MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement | 81044 | 10 bottles |

One bottle is sufficient to prepare 500 ml of medium

TABLE OF SYMBOLS

| | | | | |
|-----------------------------|---|--|--|--|
| LOT Batch code |  Do not reuse |  Manufacturer |  Contains sufficient for <n> tests |  Temperature limitation |
| REF Catalogue number |  Fragile, handle with care |  Use by |  Caution, consult accompanying documents | |

LIOFILCHEM s.r.l.

Via Scozia Zona Ind.le - 64026 Roseto D.A. (TE) - Italy

Tel. +390858930745

Fax +390858930330

Website: www.liofilchem.net

E-Mail: liofilchem@liofilchem.net

Rev.0/11.06.2006

MODIFIED LISTERIA SELECTIVE ENRICHMENT Supplement

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