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BIBLIOGRAFIA

- Van Netten, P. et al., (1989). J. of Food Microbiol. 8:299-317.
- Lovett, J., D.W. Francis, J.M. Hunt (1988) *Listeria monocytogenes* in raw milk; detection, incidence and pathogenicity. . Food Protect. 50:188-192.
- Lovett, J., A.D.Hitchins (1989) *Listeria* isolation. FDA Bacteriological Analytical Manual. 6th Ed. SUPP. SEPT. 1987 (2nd Print):29.01
- FDA Bacteriological Analytical Manual; 8th Edition, 1998. Chapter 10.

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| | | |
|--|------------|---|
| Prodotto | REF |  |
| LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF) | 81045 | 10 flaconi |

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Rev.0 / 18.07.2007

LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)

Supplement for the selective enrichment of *Listeria spp.* in food and milk samples

DESCRIPTION

LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF) is a selective supplement made of a freeze-dried mixture of nalidixic acid, cicloheximide and acriflavine to use as supplement of the culture medium LISTERIA ENRICHMENT B. BASE (Lovett) code 610326 or 620326 for the selective enrichment of *Listeria spp.* The supplement is recommended by FDA for food analysis and is recommended by IDF/FIL for the selective enrichment of *Listeria* in milk samples.

KIT CONTENTS

Each kit contains:

- 10 vials of freeze-dried **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)**
- 1 instructions sheet

PRINCIPLE OF THE METHOD

Nalidixic acid, cicloheximide and acriflavine contribute to the final medium selectivity by inhibiting the growth of most of common bacterial species non-*Listeria spp* frequently found in food. Nalidixic acid acts against Gram-negative bacteria, cicloheximide inhibits yeasts growth, acriflavine inhibits many Gram-positive bacteria.

COMPOSITION

| LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF) | | |
|---|-----------------------|------------------------------|
| | <i>Content / vial</i> | <i>Content / l of medium</i> |
| Nalidixic Acid | 20.0 mg | 40.0 mg |
| Cicloheximide | 25.0 mg | 50.0 mg |
| Acriflavine | 7.5 mg | 15.0 mg |

PROCEDURE FOR USE

1. Reconstitute aseptically the content of one vial of **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)** with 5 ml of sterile distilled water. Shake until completely dissolved, avoiding foam formation.
2. Add aseptically the entire content of one vial (5 ml) to 500 ml of medium LISTERIA ENRICHMENT B. BASE (Lovett) code 610326 or 620326 autoclaved and cooled at 45-50°C.
3. Mix with care.
4. Distribute into sterile final containers.

TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical documentation for LISTERIA ENRICHMENT B. BASE (Lovett) code 610326 or 620326.

QUALITY CONTROL

1. Control of the appearance: freeze-dried product, yellow colour .
2. Microbiological control.

Prepare the tubes using as base the medium LISTERIA ENRICHMENT B. BASE (Lovett) code 610326 or 620326 added with **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)** (1 vial in 500 ml of medium). The tubes are inoculated with the strains indicated in the microbiological control table.

Incubation condition: 24-48 h at 36 ± 1 °C.

Microbiological control:

| | Control strains | Growth |
|-------------------------------|------------------------|---------------|
| <i>Listeria monocytogenes</i> | ATCC 19111 | Good |
| <i>Listeria monocytogenes</i> | ATCC 13932 | Good |
| <i>Escherichia coli</i> | ATCC 25922 | Inhibited |
| <i>Enterococcus faecalis</i> | ATCC 29212 | Inhibited |

PRECAUTIONS

The product **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)** is classifiable as hazardous under current legislation; it is recommended that the Safety Data Sheet be consulted on its use.

LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF) is intended for use in a professional environment and must be used in the laboratory by properly trained personnel, using approved asepsis and safety methods for handling pathogenic agents.

STORAGE

Store **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)** at 2-8 °C in its original packaging. In such conditions **LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF)** will remain valid until the expiry date indicated on the label. Do not use beyond that date. Eliminate without using if there are signs of deterioration.

REFERENCES

- Van Netten, P. et al.,(1989). J. of Food Microbiol. 8:299-317.
- Lovett, J., D.W. Francis, J.M. Hunt (1988) *Listeria monocytogenes* in raw milk; detection, incidence and pathogenicity. . Food Protect. 50;188-192.
- Lovett, J., A.D.Hitchins (1989)*Listeria* isolation. FDA Bacteriological Analytical Manual. 6th Ed. SUPP. SEPT. 1987 (2nd Print):29.01
- FDA Bacteriological Analytical Manual; 8th Edition, 1998. Chapter 10.

PRESENTATION

| Product | REF | Σ |
|--|-------|----------|
| LISTERIA FOR ENRICHMENT SUPPLEMENT (FDA/IDF) | 81045 | 10 vials |

TABLE OF SYMBOLS

| | | | | |
|-----------------------------|---|--|---|--|
| LOT Batch code |  Do not reuse |  Manufacturer |  Contains sufficient for <n> tests |  Temperature limitation |
| REF Catalogue number |  Fragile, handle with care |  Use by |  Caution, consult accompanying documents | |

LIOFILCHEM Bacteriology Products

Via Scozia Zona Ind.le - 64026 Roseto D.A. (TE) - Italy

Tel. +390858930745 Fax +390858930330 Website: www.liofilchem.net E-Mail: liofilchem@liofilchem.net

Rev.0 / 18.07.2007