Chromatic™ *Bacillus cereus* Supplement

Selective enrichment supplement for detection of *Bacillus cereus*.

DESCRIPTION

Chromatic™ *Bacillus cereus* Supplement consists of two distinct supplements used for the preparation of Chromatic™ *Bacillus cereus* (ref. 610628, 620628). The complete medium is used for the isolation, differentiation and enumeration of *B. cereus* from a wide variety of foods and environmental samples in the area of food production and handling.

KIT CONTENTS

Each kit contains:

- 10 vials of liquid Chromatic™ *Bacillus cereus* Enrichment Supplement
- 10 vials of lyophilized Chromatic™ *Bacillus cereus* Selective Supplement
- 1 instructions sheet

PRINCIPLE OF THE METHOD

Selectivity is due to the combination of polymyxin B with trimethoprim. Polymyxin B inhibits most Gram-negative organisms and some Gram-positive bacteria including some bacilli other than *B. cereus*. Trimethoprim, which blocks folic acid synthesis necessary for cellular replication, is active against many Gram-positive bacteria including *S. aureus*, *Enterococcus* spp and some non-*Bacillus* spp.

COMPOSITION

		Content / vial	Content / liter of medium
Enrichment Supplement		1.5 g	3.0 g
Selective Supplement	Polymyxin B	53,000 IU	106,000 IU
	Trimethoprim	0.005 g	0.01 g

PROCEDURE FOR USE

1. Reconstitute aseptically the content of one vial of Chromatic™ *Bacillus cereus* Selective Supplement, with 5 ml of sterile distilled water.
2. Mix to complete dissolution and add aseptically to 500 ml of Chromatic™ *Bacillus cereus* along with one vial of Chromatic™ *Bacillus cereus* Enrichment Supplement. Autoclave and cool the medium at 45-50°C before the addition of supplements.
3. Mix with care and pour into Petri dishes.

TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical sheet of the medium being prepared.

QUALITY CONTROL

1. Visual inspection: lyophilized supplement appears as a whitish button, limpid solution once reconstituted; liquid supplement is transparent, viscous.
2. Microbiological control.
Prepare the medium per label directions. Inoculate the plates with the microbial strains indicated below and incubate at 30 ± 2°C for 18-24 h.

Control strains		Growth	Colony Appearance
<i>Bacillus cereus</i>	ATCC® 11778	Good	Blue-green colonies with opaque halo
<i>Bacillus subtilis</i>	ATCC® 6633	Inhibited	---
<i>Escherichia coli</i>	ATCC® 25922	Inhibited	---

WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional only and must be used by properly trained operators.

STORAGE AND TRANSPORT CONDITIONS

2-8°C away from light, until the expiry date on the label. However, our stability studies have shown that the storage or transport at 18-25°C for 4 days, or at 35-39°C for 48 hours, do not alter in any way the performance of the product. Eliminate if signs of deterioration or contamination are evident.

REFERENCES

- ISO 7932:2004. Horizontal method for the enumeration of presumptive *Bacillus cereus*.
- ISO 16140:2003. Microbiology of food and animal feeding stuffs - Protocol for the validation of alternative
- Foodborne Pathogenic Microorganisms and Natural Toxins Handbook *Bacillus cereus* and other *Bacillus* spp. (2003) U.S. Food & Drug Administration (C.F.S.A.N)
- Handbook of Culture Media for Food Microbiology (2003) Volume 37. Chapter 4. Media for *Bacillus* spp. and related genera relevant to foods. Edited by Corry, J. E. L., Curtis, G. D. W. and Baird, R. M. Publisher - Elsevier, Amsterdam.

PRESENTATION

Product	Ref.	Content
Chromatic™ <i>Bacillus cereus</i> Supplement	81097	20 vials

One vial of Enrichment Supplement and one vial of Selective Supplement are sufficient to prepare 500 ml of medium.

TABLE OF SYMBOLS

LOT Batch code	 Do not reuse	 Manufacturer	 Contains sufficient for <n> tests	 Temperature limitation
REF Catalogue number	 Fragile, handle with care	 Use by	 Caution, consult accompanying documents	

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LIOFILCHEM® S.r.l.