# Preston Supplement

Selective supplement for detection of *Campylobacter* spp according to ISO 10272-1.

## DESCRIPTION

Preston Supplement is a lyophilized selective supplement used in combination with Horse Blood Lysed (ref. 83397) for the enrichment of *Campylobacter* from food, animal feeding and environmental samples. Nutrient Broth No 2 ISO 10272-1 (ref. 610393, 620393) is used as basic medium. The complete medium (Preston broth) is recommended in the ISO method for samples with low numbers of campylobacters and high level of background microflora (detection procedure B).

## KIT CONTENTS

Each kit contains:

- 10 vials of freeze-dried Preston Supplement;
- 1 instruction sheet.

## PRINCIPLE OF THE METHOD

Preston Supplement contains antibiotics inhibiting contaminant microflora.

## COMPOSITION

	Content / vial	Content / liter of medium
Polymyxin B Sulfate	2 500 IU	5 000 IU
Rifampicin	5.0 mg	10.0 mg
Trimethoprim Lactate Salt	5.0 mg	10.0 mg
Amphotericin B	5.0 mg	10.0 mg

## PROCEDURE FOR USE

1. Aseptically reconstitute the content of one vial of Preston Supplement with 5 ml of ethanol 95%, mix to complete dissolution
2. To Nutrient Broth No 2 (470 ml), autoclaved and cooled to 45-50°C, add the blood (25 ml), then the antibiotic solution ( 5ml) contained in the vial.
3. Mix well and dispense into final containers.

## TECHNIQUE AND INTERPRETATION OF THE RESULTS

Refer to the technical sheet of the medium being prepared.

## QUALITY CONTROL

1. Control of the appearance: freeze-dried orange in colour.
2. Microbiological control.  
Prepare the medium per label directions. Inoculate the medium with the strains indicated below and incubate at  $41.5 \pm 1^\circ\text{C}$  for  $24 \pm 2$  h under microaerobic atmosphere.

### Control strains

### Specification

<i>Campylobacter jejuni</i> + <i>Escherichia coli</i>	WDCM 00156 WDCM 00012	>10 characteristic colonies on mCCD agar
<i>Escherichia coli</i>	WDCM 00012	Totally inhibited on TSA

## WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

## STORAGE

2-8°C in its original packaging. Keep away from sources of heat and avoid excessive changes of temperature. Use until the expiry date indicated on the label. Eliminate without using if there are signs of deterioration.

## REFERENCES

- ISO 10272-1:2017. Microbiology of the food chain – Horizontal method for detection and enumeration of *Campylobacter* spp. – Part 1: Detection method.

## PRESENTATION

Product	Ref.	Contents
Preston Supplement ISO 10272-1	81100	10 vials

One vial is sufficient to prepare 500 ml of medium.

## TABLE OF SYMBOLS

 Batch code	 Do not reuse	 Manufacturer	 Contains sufficient for <n> tests	 Temperature limitation
 Catalogue number	 Fragile, handle with care	 Use by	 Caution, consult accompanying documents	

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**LIOFILCHEM® S.r.l.**