## Easy Dry™ TBX

Chromogenic medium for detection and enumeration of *E. coli*, according to ISO 16649-1.

### DESCRIPTION

Liofilchem Easy Dry™ are absorbent pads impregnated with a sterile, dehydrated culture medium. Each pad is preplated in a Petri dish and is immediately ready to use after pouring sterile distilled or deionized water on it. Easy Dry™ are optimal for the examination of large sample volumes by the membrane filter method.

Easy Dry™ TBX is a chromogenic selective medium used for the isolation, identification and enumeration of  $\beta$ -glucuronidase-positive *Escherichia coli* in foodstuffs, animal feed and water.

This medium complies with the recommendations of ISO 16649-1.

| TYPICAL FORMULA            | (g/l) |
|----------------------------|-------|
| Enzymatic Digest of Casein | 20.0  |
| Bile Salts No. 3           | 1.5   |
| X-Glucuronide              | 0.075 |
| Final pH 7.2 ± 0.2 at 25°C |       |

### METHOD PRINCIPLE

Enzymatic digest of casein provides amino acids, nitrogen, carbon, vitamins and minerals for organisms growth. Bile salts act as selective agent inhibiting most Gram-positive bacteria. 5-bromo-4-chloro-3-indolyl- $\beta$ -D-glucuronide (BCIG) is the chromogenic substrate cleaved by the enzyme  $\beta$ -glucuronidase of *E. coli*.

### PREPARATION

1. Cut open a bag and remove the number of Easy Dry™ plates needed.
2. Moisten the pad contained in the Petri dish with 2.2 ml of sterile distilled or deionized water.
3. Wait 5 minutes before using.

### TEST PROCEDURE

For the recovery of sub-lethally injured *E. coli*, a resuscitation step is recommended by ISO 16649-1:

Place a membrane filter onto the dried surface of a plate of Mineral-Modified Glutamate Agar (MMGA). Spread 1 ml of the test sample evenly over the whole membrane surface and incubate at 37°C for 4 ± 1 h.

Under the same conditions, using decimal dilutions of the test sample or of the initial suspension, two plates per dilution are inoculated.

After resuscitation, transfer the membrane from MMGA to a plate containing a just rehydrated pad of Easy Dry™ TBX. Incubate aerobically at 44°C for further 18-24 h.

### INTERPRETING RESULTS

After incubation observe the color of the colonies and interpret the results as indicated in the ID table.

#### ID Table.

| Microorganism   | Typical colony color |
|---|----------------------|
| $\beta$ -glucuronidase-positive <i>Escherichia coli</i>     | Blue to blue-green   |
| $\beta$ -glucuronidase-negative bacteria (if not inhibited) | White to green-beige |

Count the typical colony-forming units (CFU) of  $\beta$ -glucuronidase-positive *Escherichia coli* in each plate containing less than 150 typical CFU and less than 300 total (typical and non-typical) CFU.

Note:  $\beta$ -glucuronidase-negative *E. coli* strains form colorless colonies on this medium, e.g. *E. coli* O157, or they cannot grow at elevated temperature of 44°C, e.g. *E. coli* O157:H7.

Typical reactions for confirmation of *E. coli* are a positive indole-reaction as well as a negative oxidase- and catalase-reaction.

### APPEARANCE OF THE MEDIUM

Whitish pad. Yellowish once rehydrated.

**STORAGE**

Store at 10-25°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

**SHELF LIFE**

2 years.

**QUALITY CONTROL**

Plates are inoculated with the microbial strains indicated in the QC table.

Inoculum for productivity: 50-100 CFU.

Inoculum for selectivity: 10<sup>4</sup>-10<sup>6</sup> CFU.

Incubation conditions: aerobically at 44 ± 1°C for 18-24 hours.

**QC Table.**

| Microorganism                 |            | Growth    | Specification                 |
|-------------------------------|------------|-----------|-------------------------------|
| <i>Escherichia coli</i>       | WDCM 00013 | Good      | Blue to blue-green colonies   |
| <i>Escherichia coli</i>       | WDCM 00202 | Good      | Blue to blue-green colonies   |
| <i>Enterococcus faecalis</i>  | WDCM 00009 | Inhibited | ---                           |
| <i>Citrobacter freundii</i>   | WDCM 00006 | Good      | White to green-beige colonies |
| <i>Pseudomonas aeruginosa</i> | WDCM 00025 | Good      | White to green-beige colonies |

**WARNING AND PRECAUTIONS**

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

**DISPOSAL OF WASTE**

Disposal of waste must be carried out according to national and local regulations in force.

**BIBLIOGRAPHY**

- EN ISO 11133:2014. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media.
- ISO 16649:2001. Microbiology of food and animal feeding stuffs – Horizontal method for the enumeration of β-glucuronidase-positive *Escherichia coli* – Part 1: Colony-count technique at 44°C using membranes and 5-bromo-4-chloro-3-indolyl-β-D-glucuronide.
- Ogden I.D., and A.J. Watt (1991) An evaluation of fluorogenic and chromogenic assays for the direct enumeration of *E. coli*. Letters in App. Microbiol. 13:212-215.
- Delisle G.L., and A. Ley (1989) Rapid detection of *E. coli* in urine samples by a new chromogenic β-Dglucuronidase assay. J. Clin. Microbiol. 27:778-779.

| PRESENTATION  | Contents                  | Ref.             |
|---------------|---------------------------|------------------|
| Easy Dry™ TBX | 60 mm ready-to-use plates | 100 plates 87503 |

**TABLE OF SYMBOLS**

|                             |   |   |   |   |
|-----------------------------|---|---|---|---|
| <b>LOT</b> Batch code       |  Keep away from sunlight |  Manufacturer                      |  Use by                               |  Fragile, handle with care |
| <b>REF</b> Catalogue number |  Temperature limitation  |  Contains sufficient for <n> tests |  Caution, consult Instruction For Use |  Do not reuse              |

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