Easy Dry™ Pseudomonas CN

Selective medium for detection and enumeration of *Pseudomonas aeruginosa* in water, according to ISO 16266.

DESCRIPTION

Liofilchem® Easy Dry™ are sterile, dehydrated culture media, impregnated upon absorbent pads. Each pad is preplated in a Petri dish and is immediately ready-to-use after being moistened with sterile distilled or deionized water. Easy Dry™ are optimal for the examination of large sample volumes by the membrane filter method.

Easy Dry™ Pseudomonas CN is a selective isolation medium used for the detection and enumeration of *Pseudomonas aeruginosa* in water, based on the promotion of pigment production, following the procedure given by EN ISO 16266.

TYPICAL FORMULA

Gelatin Peptone

Casein Hydrolysate

Potassium Sulfate

Magnesium Chloride

Glycerol

Cetrimide

Nalidixic Acid

Final pH 7.1 ± 0.2 at 25°C

Formula is based on the corresponding agar medium, adjusted to meet specific performance requirements.

METHOD PRINCIPLE

Gelatin peptone and casein hydrolysate provide amino acids, nitrogen, carbon, vitamins and minerals.

Potassium sulfate and magnesium chloride stimulate pyocyanin production. Glycerol is a carbon and energy source. Cetrimide (hexadecyltrimethyl ammonium bromide) is a quaternary ammonium compound that inhibits a wide variety of organisms. Nalidixic acid is effective against both Gram-positive and Gram-negative bacteria other than *Pseudomonas aeruginosa*.

PREPARATION

1. Cut open a bag and remove the number of Easy Dry™ plates needed.
2. Moisten the pad contained in the Petri dish with 2.2 ml of sterile distilled or deionized water.
3. Wait 5 minutes before using.

TEST PROCEDURE

Filter the sample through a filter membrane (0.45 µm pore diameter). Transfer the membrane onto a plate containing a just rehydrated pad. Incubate aerobically at 36 ± 2°C for 40-48 hours.

INTERPRETING RESULTS

Examine the membrane for growth and fluorescence under UV light at 360 nm after 20-24 h and 40-48 h.

- Count all colonies producing the green-blue pigment as confirmed *Pseudomonas aeruginosa*.
- Non-pyocyanin producing colonies that fluoresce under UV should be confirmed by the acetamide test.
- All other reddish-brown non-pigmented colonies that do not fluoresce should be submitted to further testing such as ammonia production from acetamide, oxidase production and fluorescence on King's B medium.

For confirmation, select as many colonies as possible (all colonies if possible). Streak each colony onto a Nutrient Agar (NA) plate and incubate at 36 ± 2°C for 22 ± 2 h. Select one well isolated colony from each NA plate for the confirmatory tests.

APPEARANCE OF THE MEDIUM

Whitish pad. Yellowish once rehydrated.

STORAGE

Store at 10-25°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

SHELF LIFE

2 years.

QUALITY CONTROL

The medium is inoculated with the microbial strains indicated in the QC table.

Inoculum for productivity: 50-100 CFU.

Inoculum for selectivity: 10⁴-10⁶ CFU.

Incubation conditions: aerobically at 36 ± 2°C for 40-48 hours.

QC Table.

Microorganism	Growth
<i>Pseudomonas aeruginosa</i>	WDCM 00024
<i>Pseudomonas aeruginosa</i>	WDCM 00025
<i>Enterococcus faecalis</i>	WDCM 00009
<i>Escherichia coli</i>	WDCM 00012

WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to national and local regulations in force.

BIBLIOGRAPHY

See the references at the end of this document.

PRESENTATION	Packaging	Ref.
Easy Dry™ Pseudomonas CN	100 pads	87505

TABLE OF SYMBOLS

See the table of symbols at the end of this document.

This document is available from the online Support Center:

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liofilchem.com/ifu-sds

BIBLIOGRAPHY / BIBLIOGRAFÍA

1. EN ISO 11133:2014+Amd1:2018. Microbiology of food, animal feed and water – Preparation, production, storage and performance testing of culture media.
2. EN ISO 16266:2008. Water quality – Detection and enumeration of *Pseudomonas aeruginosa* by membrane filtration.
3. Goto S. and S. Enomoto (1970) Nalidixic acid cetrimide agar. A new selective plating medium for the selective isolation of *P. aeruginosa*. Jpn. J. Microbiol. 14:65.
4. Brown V.L. and E.J.L. Lowbury (1965) Use of an improved Cetrimide Agar Medium and of culture methods for *P. aeruginosa*. J. Clin. Pathol. 18:752.
5. King E.O., M.K. Ward and D.E. Raney (1954) Two simple media for the demonstration of pyocyanin and fluorescin. J. Lab. Clin. 44, 301.

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