Easy Dry™ MacConkey

Selective and differential medium for detection of Enterobacteriaceae in water, food and other materials, according to USP/EP/JP.

DESCRIPTION

Liofilchem Easy Dry™ are absorbent pads impregnated with a sterile, dehydrated culture medium. Each pad is preplated in a Petri dish and is immediately ready to use after pouring sterile distilled or deionized water on it. Easy Dry™ are optimal for the examination of large sample volumes by the membrane filter method.

Easy Dry™ MacConkey is a slightly selective medium giving excellent differentiation between lactose-fermenting and lactose-nonfermenting Gram-negative enteric bacilli, from foodstuffs, waste water, pharmaceutical, and industrial sources.

This medium conforms to harmonized USP/EP/JP requirements for the detection of *E. coli* in nonsterile pharmaceutical products.

TYPICAL FORMULA	(g/l)
Pancreatic Digest of Gelatin	17.0
Peptone from Meat	1.5
Peptone from Casein	1.5
Lactose	10.0
Sodium Chloride	5.0
Bile Salts	1.5
Neutral Red	0.03
Crystal Violet	0.001
Final pH 7.1 ± 0.2 at 25°C	

METHOD PRINCIPLE

Pancreatic digest of gelatin and peptones from meat and casein provide amino acids, nitrogen, carbon, vitamins and minerals for organisms growth. Lactose is the fermentable carbohydrate. Sodium Chloride maintains the osmotic balance of the medium. Bile salts and crystal violet are the selective agents, inhibiting Gram-positive organisms and allowing Gram-negative bacteria to grow. Neutral red is the pH indicator.

PREPARATION

1. Cut open a bag and remove the number of Easy Dry™ plates needed.
2. Moisten the pad contained in the Petri dish with 2.2 ml of sterile distilled or deionized water.
3. Wait 5 minutes before using.

TEST PROCEDURE

Filter the sample trough a filter membrane (0.45 µm pore diameter). Transfer the membrane onto a plate containing a just rehydrated pad. Incubate aerobically at 30-35°C for 18-72 h.

INTERPRETING RESULTS

Lactose-nonfermenting organisms, such as *Salmonella*, *Shigella* and *Proteus* spp, form colorless or clear colonies. Lactose-fermenting organisms, such as *E. coli* and *Klebsiella* spp, grow as pink to red colonies with or without a zone of precipitated bile. Enterococci, Staphylococci and other Gram-positive bacteria are partially or completely inhibited.

To confirm the presence of *E. coli*, transfer the membrane and the colonies onto a pad of Easy Dry™ TBX (ref. 87503), just rehydrated and pre-warmed to 44°C, for the detection of β-glucuronidase activity. More than 94 % of *E. coli* strains are β-glucuronidase-positive and will grow as blue to blue-green colonies on TBX Agar. Typical reactions for detection of *E. coli* are a positive indole-reaction as well as a negative oxidase- and catalase-reaction.

APPEARANCE OF THE MEDIUM

Pinkish pad. Slightly darker once rehydrated.

STORAGE

Store at 10-25°C away from light. Do not use the product beyond its expiry date on the label or if product shows any evidence of contamination or any sign of deterioration.

SHELF LIFE

2 years.

QUALITY CONTROL

Plates are inoculated with the microbial strains indicated in the QC table.

Inoculum for productivity: 50-100 CFU.

Inoculum for selectivity: 10⁴-10⁶ CFU.

Incubation conditions: aerobically at 30-35°C for 18-72 hours.

QC Table.

Microorganism		Growth	Specification
<i>Salmonella Typhimurium</i>	ATCC® 14028	Good	Colorless colonies
<i>Shigella flexneri</i>	ATCC® 12022	Good	Colorless colonies
<i>Proteus mirabilis</i>	ATCC® 12453	Good	Colorless colonies
<i>Escherichia coli</i>	ATCC® 8739	Good	Pink colonies with bile ppt
<i>Klebsiella pneumoniae</i>	ATCC® 13883	Good	Pink colonies
<i>Enterococcus faecalis</i>	ATCC® 29212	Partially to completely inhibited	Very small, opaque colonies
<i>Staphylococcus aureus</i>	ATCC® 25923	Inhibited	---

WARNING AND PRECAUTIONS

The product does not contain hazardous substances in concentrations exceeding the limits set by current legislation and therefore is not classified as dangerous. It is nevertheless recommended to consult the safety data sheet for its correct use. The product is intended for professional use only and must be used by properly trained operators.

DISPOSAL OF WASTE

Disposal of waste must be carried out according to national and local regulations in force.

BIBLIOGRAPHY

- European Pharmacopoeia 6.5 (2009). 2.6.13 Microbiological examination of non-sterile products: Test for specified microorganisms.
- United States Pharmacopoeia 32 NF 27 (2009). <62> Microbiological examination of non-sterile products: Test for specified microorganisms.
- Japanese Pharmacopoeia 4.05 (2008). Microbiological examination of non-sterile products: Test for specified microorganisms.
- Murray, Baron, Jorgensen, Landry and Pfaller ed. (2007) Manual of clinical microbiology, 9th ed. American Society for Microbiology, Washington, D.C.
- MacConkey A. (1905) Lactose-fermenting bacteria in faeces. J. Hygiene 8:333-379.

PRESENTATION	Contents	Ref.
Easy Dry™ MacConeky	60 mm ready-to-use plates	100 plates

TABLE OF SYMBOLS

LOT Batch code	 Keep away from sunlight	 Manufacturer	 Use by	 Fragile, handle with care
REF Catalogue number	 Temperature limitation	 Contains sufficient for <n> tests	 Caution, consult Instruction For Use	 Do not reuse

**LIOFILCHEM® s.r.l.**Via Scozia zona ind.le, 64026 Roseto degli Abruzzi (Te) Italy
Tel. +39 0858930745 Fax +39 0858930330 www.liofilchem.net liofilchem@liofilchem.net

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