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Fare riferimento a Standard Methods for the Examination of Dairy Products.

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E-Mail: liofilchem@liofilchem.net

Rev.4 / 22.12.2017

Suspension of *Bacillus subtilis* (ATCC 6633) spores
for determination of penicillin in milk, food and biological fluids.

DESCRIPTION

SUSPENSION SPORE BAS E5/E6/E7/E8/E9/E10 is a standard and stabilized suspension of *Bacillus subtilis* ATCC 6633 used to detect the presence of penicillin up to 0.05 UI in milk and of other antibiotics according to Standard Methods for the Examination of Dairy Products.

CONTENT OF THE PACKAGE

Each package contains:

- Suspension of *Bacillus subtilis* ATCC 6633 spores
- Instruction sheet

COMPOSITION

SUSPENSION SPORE BAS E5/E6/E7/E8/E9/E10 contains spores of *Bacillus subtilis* ATCC 6633 in distilled water in predefined concentrations: E5=1-5x10⁵CFU/ml; E6=1-5x10⁶CFU/ml; E7=1-5x10⁷CFU/ml; E8=1-5x10⁸CFU/ml; E9=1-5x10⁹CFU/ml; E10=1-5x10¹⁰CFU/ml.

TECHNIQUE AND RESULTS INTERPRETATION

Refer to Standard Methods for the Examination of Dairy Products.

QUALITY CONTROL

1. Appearance: turbid solution, milky.
2. Sensitivity to penicillin:

Antibiotic concentration	Inhibition halo
PENICILLIN G 217,1 IU	38-32 mm
PENICILLIN G 87 IU	30-36 mm
PENICILLIN G 0,86 IU	21-27 mm
PENICILLIN G 0,08 IU	13-19 mm
PENICILLIN G 0,052 IU	11-17 mm

PRECAUTIONS

The product **SUSPENSION SPORE BAS E5/E6/E7/E8/E9/E10** is not classified as hazardous by current legislation. The product is designed for a professional environment and must be used only by properly trained operators.

STORAGE

Store the product at 2-8°C: in these conditions it maintains its validity until the expiry date indicated on the label.

BIBLIOGRAFIA

1. *Standard Methods for the Examination of Dairy Products*, APHA, Washington, D.C., 14th Ed. 1978.

PRESENTATION

PRODUCT	REF	PACKAGING
SUSPENSION SPORE BAS E5	82015	1 vial x 10 mL
SUSPENSION SPORE BAS E6	82014	1 vial x 10 mL
SUSPENSION SPORE BAS E7	82013	1 vial x 10 mL
SUSPENSION SPORE BAS E8	82012	1 vial x 10 mL
SUSPENSION SPORE BAS E8	82212	1 vial x 2 mL
SUSPENSION SPORE BAS E8	82112	1 bottle x 100 mL
SUSPENSION SPORE BAS E9	82020	1 vial x 10 mL
SUSPENSION SPORE BAS E10	82019	1 vial x 10 mL

TABLE OF SYMBOLS

 Manufacturer	 Contains sufficient for <n> tests	 Temperature limitation
 Catalogue number	 Fragile, handle with care	 Caution, consult accompanying documents
 Use by	 Batch code	 Do not reuse

LIOFILCHEM® S.r.l. Microbiology Products

Via Scozia Zona Ind.le - 64026 Roseto D.A. (TE) - Italy

Tel. +390858930745

Fax +390858930330

Website: www.liofilchem.net

E-Mail: liofilchem@liofilchem.net

Rev.4 / 22.12.2017